


Programming for Modes for BX-Series Scales

1. Turn the scale ON and once the display shows the ROM version number, press the  key.




*Note: The version number may vary.

2. The display will then show:




3. Press , , , , and then .

4. The display should look like this:

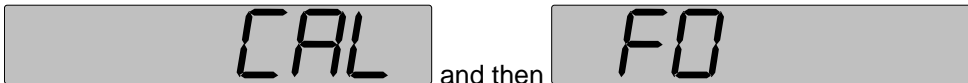

5. Press the  key repeatedly until the display reads F17:



6. Now press the  key and the display will look like this:




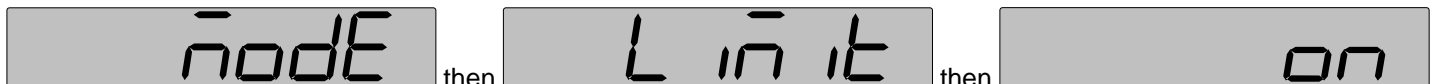
7. You can use the  key to enable or disable the Weigh1 Mode (1/3000; 1/6000; or 1/12,000 Resolution.)


8. Now press the  key and the display will look like this:



9. You can use the  key to enable or disable the Weigh2 Mode (1/30,000; 1/60,000; or 1/120,000 Resolution.)


10. Now press the  key and the display will look like this:



11. You can use the  key to enable or disable the Limit Mode.


12. Now press the  key and the display will look like this:



13. You can use the  key to enable or disable the Hold Mode.

14. Now press the  key and the display will look like this:



15. Press  the key and the scale will do the Display Self-Test and then go to zero.

If any problems persist, please contact the Atron Systems Service Department at 973-227-8882.