
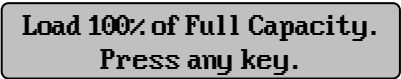



LS-100 Quick Calibrate Instructions, kg

1. Make sure that the scale is plugged into a proper outlet and that the *Power Switch* is **OFF**.
2. Locate the *CAL Switch* which is under the platter on the front left side of the scale.
You may need to remove the *CAL Plate* to expose the *CAL Switch*.
3. Press and hold down the *CAL Switch* as you place the *Power Switch* in the **ON** position.
4. Once the displays read , release the *CAL Switch*.
5. Make sure that there is nothing on the platter. Press , and .
6. Once the LCD display reads ,
7. You can press the key to toggle the percentage between 100%, 50%, 20%, and 10%.
8. Place the correct load on the platter based on the chart below:

Percentage	Calibration Weight
100 %	30 kg (or 66.14 lb)
50 %	15 kg (or 33.07 lb)
20 %	6 kg (or 13.23 lb)
10 %	3 kg (or 6.61 lb)

9. Once the weight is stable press .
10. After about 5 seconds the scale will beep again and then the LCD display will show:

Calibrated!

11. Remove all weight from the platter and then press .
12. Press , , and then .
13. The scale displays should show .
14. You are done.

If further problems persist, please contact the Atron Systems Service Department at 973-227-8882.