


PC-100 Successive Tare Programming

1. Make sure that the scale is plugged into a proper outlet, the platter is properly seated, there is nothing on the platter, and that the *Power Switch* is **OFF**.
2. Locate the *CAL Switch* which is under the platter on the front left side. If necessary, remove the *CAL Plate* to expose the *CAL Switch*.

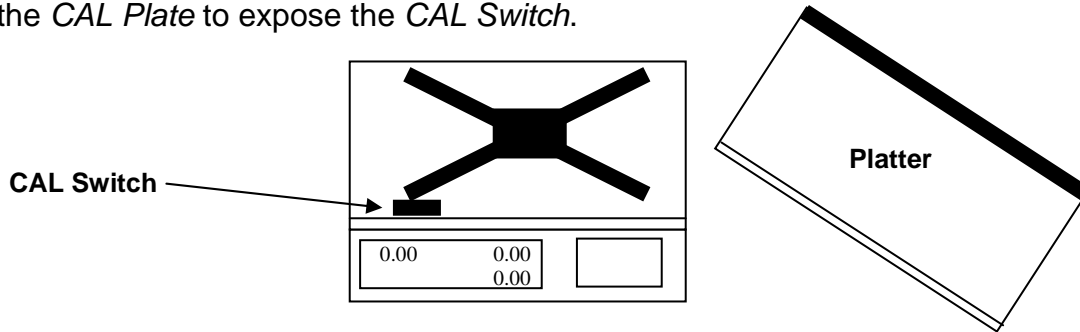

3. Press and hold down the *CAL Switch* as you place the *Power Switch* in the **ON** position.


4. The displays will read: *SPAn SEt*.

5. Press **TARE** **TARE** and then the display will read: *Code*.


6. Now press **C** and the displays will now read: *Code 000*.

7. Press **7** **7** **5** **5** **9** and the displays will now read: *Code 77559*.

8. Press the ***** key 2 times. The displays will read: *PR-R SEt 1* and the *Price Display* will have a hexadecimal number in it. Now you need to enter the following key strokes:

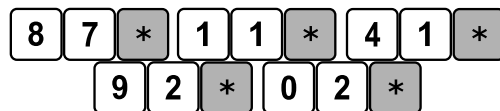

Now the displays will read *End* and then read *SPAn SEt*.

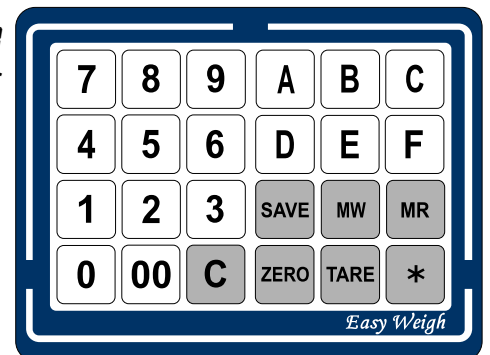

9. Go to step 4 of ***PC-100 Quick Calibrate Instructions*** and complete those instructions using full capacity weights (60 pounds.) If you encounter problems with the ***PC-100 Quick Calibrate Instructions***, it may be necessary to follow the ***PC-100 Span & Gain Calibration Instructions***.

For kilogram setup & decimal place change, see ***PC-100 kg/lb Unit Programming*** instructions.
If further problems persist, please contact the Atron Systems Service Department at 973-227-8882.